Vidrio Industrial POBEL, S.A.


Bomba de vacío
Vacuum pump

BOVT 0350

MANUAL DE INSTRUCCIONES
INSTRUCTIONS MANUAL

C/ Luis I, s/n – Almacén 12 – Pol. Ind. Vallecas IV
28031 Madrid (ESPAÑA)
Tel. (34) 91 380 33 18 – Fax (34) 91 380 32 16
E-mail: pobel@pobel.com – Web: www.pobel.com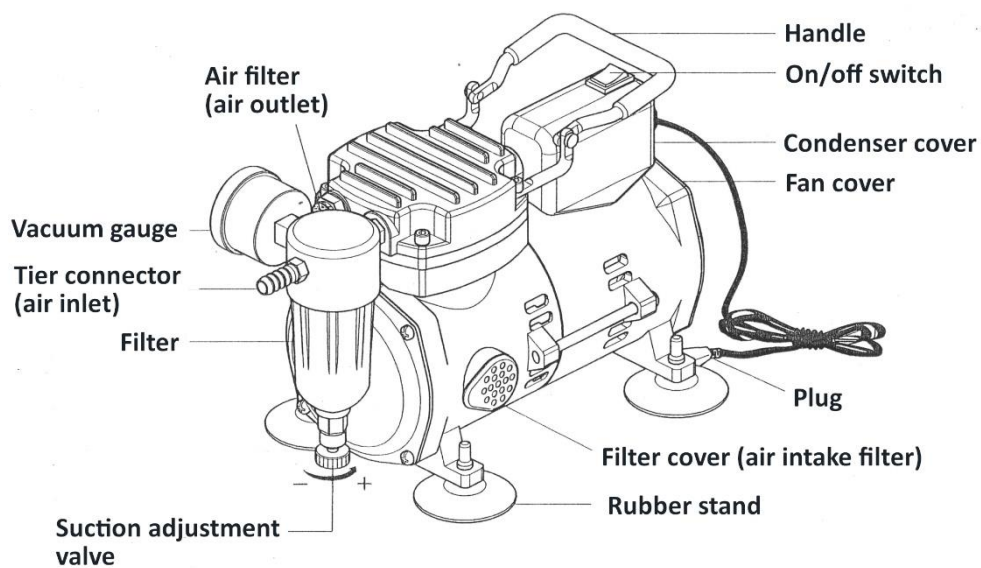

PRECAUTIONS

1. Please check the voltage and spec. label on the machine.
2. Carefully unpack the machine and check if any part is damaged during transportation. Any damage should be reported immediately to the service agent.
3. Plug the vacuum pump into mains supply, making sure the voltage is correct before you turn on the machine.
4. Work in a well ventilated, clean area and away from wet surfaces.
5. The vacuum pump has a thermal switch that automatically shuts off when overheated or lacking voltage.
6. Adjust suction by turning the suction adjustment valve clockwise for less suction, counter-clockwise for more suction.
7. Do not touch the vacuum pump during or soon after use as it may become hot.
8. Do not use any lubricants.
9. For best performance, the machine should not run continuously over one hour and should be allowed to cool down.
10. Improper operation and dismantling of the parts might damage the vacuum pump. Damaged machine must be sent to the service agent for replacement or repair to avoid hazard. Do not attempt to fix the machine without consulting the service agent first.
11. Do not use for suction of fluids or gases with high moisture content.


PROBLEM SHOOTING

Machine cannot operate

1. Check if plug is inserted correctly onto the receptacle.
2. Check if the occurrence is due to overheat, which leads to automatic shut off. Cool off for 30 minutes before turning the power back on.
3. Check if the On/off switch is turned on.

Suction reduced

1. Check for air leakage: check if all connections are sealed tight.
2. Check if the suction adjustment valve is closed tight.
3. Check if unintended objects are suctioned into the unit, causing blockage.
4. Check if voltage applied is too low.

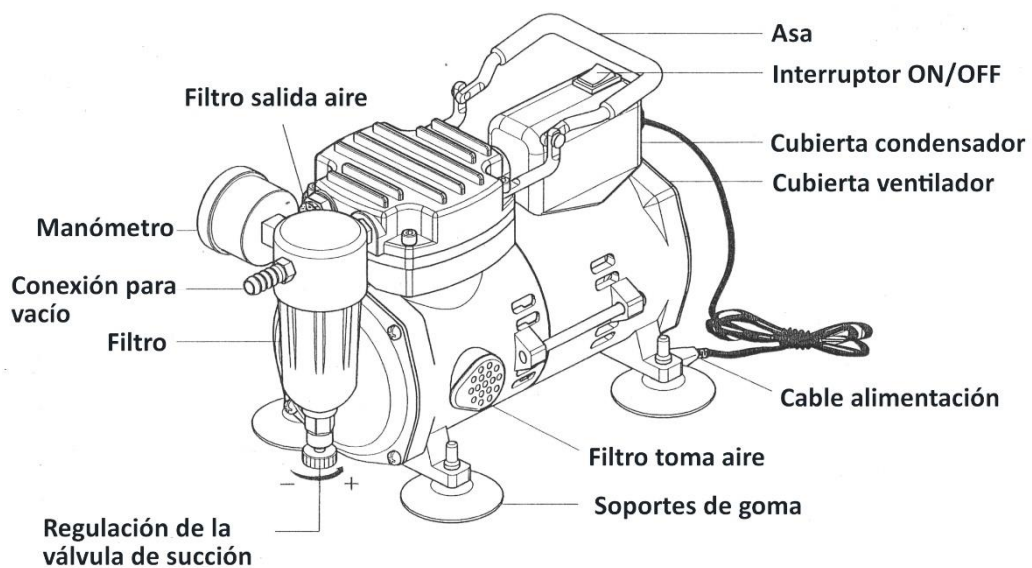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