**Vidrio Industrial POBEL, S.A.**

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## 1.1.GENERAL INFORMATION

- The package must be handled carefully. When you unpackage it, you should check that what it contains inside is what is indicated in the Packing List section. If any component is missing or otherwise damaged, contact the dealer immediately.
- Please, read the instructions before installing the apparatus. The instruction manual will be an inseparable part of the product; For this reason, keep it in good condition for future reference.
- The modification of the components of the product transgresses the directive of use 2009/104 / CE; and therefore: the manufacturer is not responsible for the damages that this brings as damage.
- The product must be installed on a horizontal surface and never in places where there are vapors of contaminating or flammable substances.
- Please, dispose of at least the first 4 liters of distilled water before using the equipment. • Conservative packaging during the warranty period.

**¡ATTENTION! IT WILL NOT BE ADMITTED TO REPAIR ANY DISTILLER THAT IS NOT PROPERLY CLEAN.**

### **IMPORTANT:**

**RINSE WITH PLENTY OF WATER TO REMOVE SALT FROM INSIDE THE CONDENSER.**

## 1.2.PACKING LIST

Description	Quantity
Boiler	1
Condenser	1
Level	1
Metallic support base	1
Water feed hose	1
Filter	1
Set of silicone tubes	1
Instructions manual	1

## 1.3.TECHNICAL SPECIFICATIONS

Reference	Length (mm)	Width (mm)	Depth (mm)	Weight (Kg)
0100710710	510	590	170	5,2

Capacity (L/H)	Water consumption (L/MIN)	Power (W)	Voltage (V)	Electric current (A)
4	1	2800	230	12.2

## 1.4.INSTALLATION

According to the photos, make the installation following these steps:

1. Join the level with the boiling body using the GL -25 thread.

COMMUNICATION BETWEEN LEVEL AND BOILER

2. Connect the gray hose to the water supply of the equipment. The other end of the hose, with the filter previously installed, will be connected to a  $\frac{3}{4}$ " tap..

3. Make sure that the seals (JV1 and JV2) are placed in the vertical tube, introduce the refrigerant vertically downwards.

JV1

JV2

4. Screw the GL-18 nut connecting the level inlet with the upper outlet or coolant outlet.

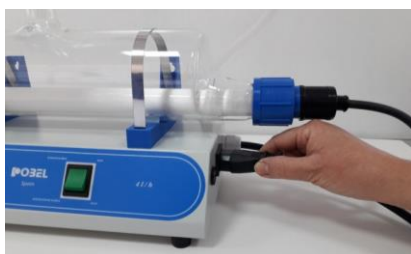
5. The pre-installed nut GL-18 on the support base joins by screwing into the cooling water inlet.


6. Carefully insert the heating element into the boiler and with the N42 nut screw it firmly until it stops.


7. Join the heating element with the distiller by inserting the pin into the female wall base (pin) of the metal base.

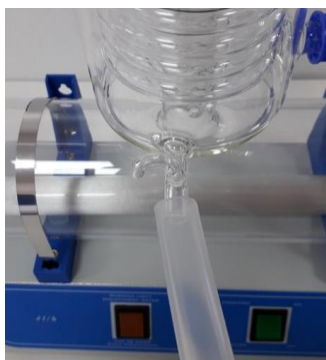8. Connect the distilled water outlet tube to the lower and front condenser inlet.

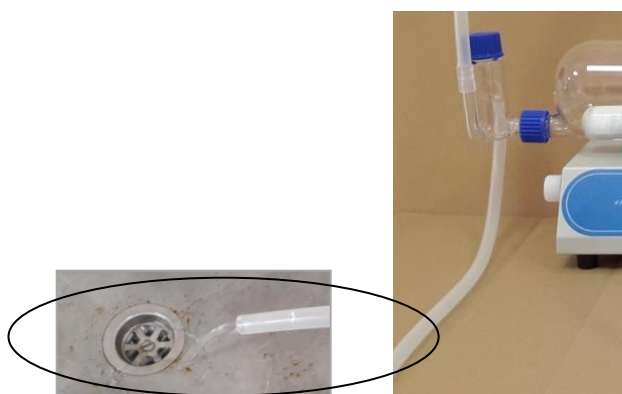
9 The level outlet pipe should be placed in a position such that, the excess of water falls freely outside in a sink.

10. Finally, connect the device to the power supply through the connection cable.

## 1.5.START-UP
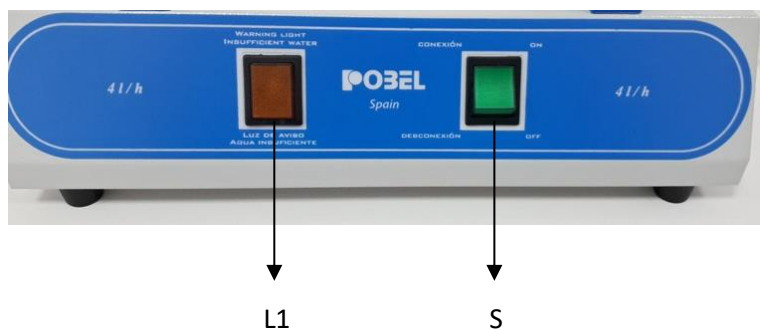
### FRONT PANEL

### IMPORTANT:

BEFORE STARTING THE COMMISSIONING VERIFY THAT THE INSTALLATION HAS BEEN MADE CORRECTLY.

In order to start the distillation, please, follow these steps:

Open the water supply tap by allowing more than 60 L/H or 1L/MIN; until the water completely covers the heating resistance.

Press the main switch (S) located on the front panel; The green light will illuminate. In the event that the yellow light (L1) lights up, we will have an insufficient water supply. If this occurs, the Flow of water must be increased until the yellow light turns out.

After approximately seven minutes, the boiling process will begin and the heating resistance acquires a visible reddish color. The distilled water that is obtained will flow through the tube intended for it. See step 8 of the installation section.

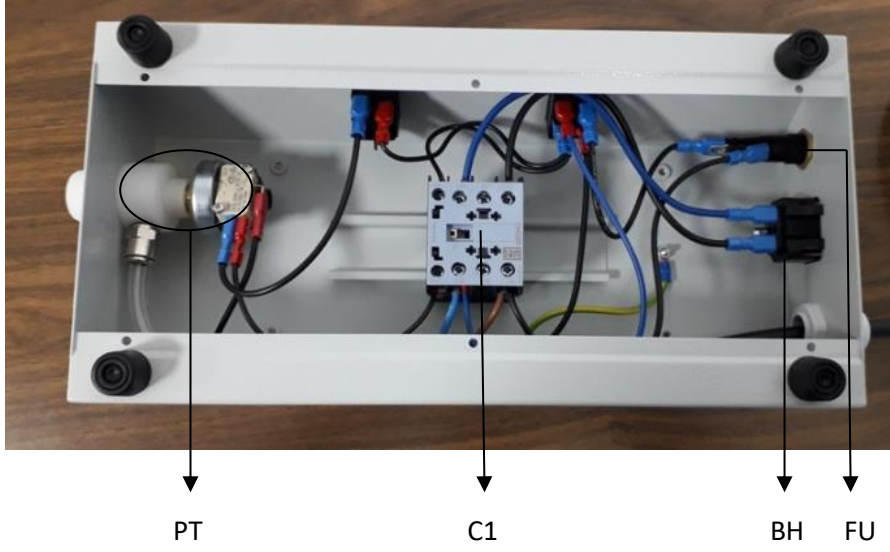
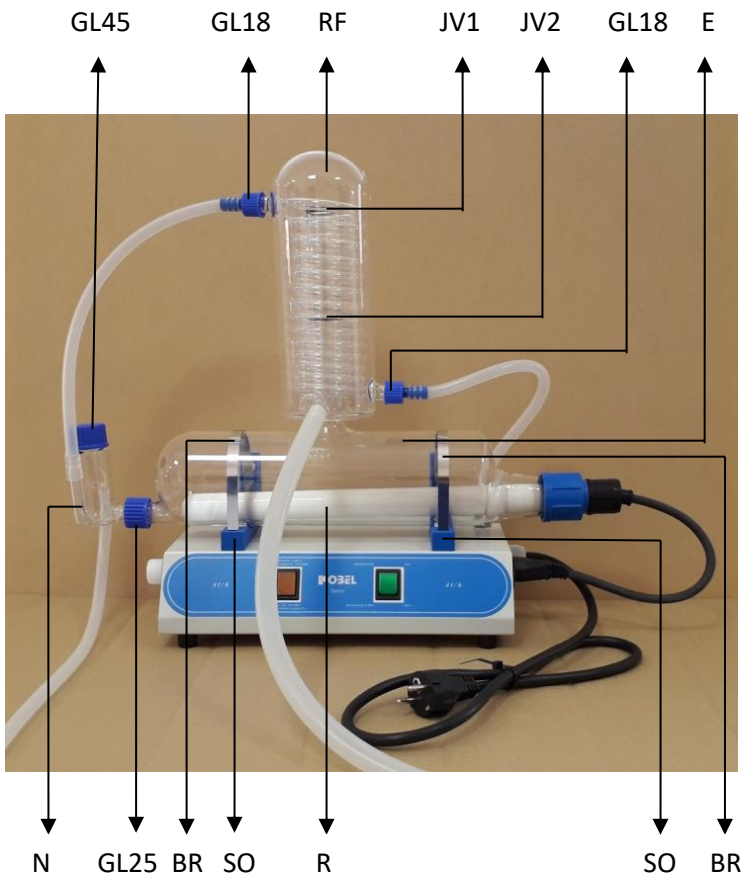
During the distillation, check that the drainage is always carried out as described in step 9 of the installation section, since right drainage of the excess of water free of obstacles guarantees to maintain the adequate level of water inside the boiler.

## **1.6.MAINTENANCE**

**ATTENTION! ALWAYS DISCONNECT THE POWER SUPPLY BEFORE CARRYING OUT THE CLEANING OPERATION.**

With the use, the lime may be deposited in the distiller, a film of this substance being formed that will decrease the heating resistance performance. Therefore, the deposited lime must be cleaned with 20% hydrochloric acid or with a commercial anti-lime product; which is poured through the inlet of the level; leaving a few minutes before rinsing properly with enough water. The frequency of the cleaning depends directly on the hardness of the water being used in the distillation process.

1.7.SPARE PARTS


DESCRIPTION	Fig.	CAT. NUMBER	QUANTITY
Quartz heater	R	0100721548	1
Flange + support f/water still	BR	0100702726	2
Blue plastic base	SO	YP.DSOPC00700	2
Viton® o-ring	JV1	YP.DJUVI21035	1
Viton® o-ring	JV2	YP.DJUVI27036	1
Glass level	N	0100919936	1
Boiler	E	0100710708	1
Condenser	RF	0100701700	1
Nut	GL25	0200161836	1
Cap	GL45	0200162060	1
Nut	GL18	0100100603	2
Silicone washer f/GL18	--	YP.CJUSI07142	2
Silicone washer f /GL25	--	YP.DJUPL24125	1
Straight plastic hose f/GL18	--	YP.DOLIV00001	2
Circuit breaker	FU	YE.DINTE00016	1
Socket	BH	YE.DCLAH00002	1
Mini contactor	C1	YE.DCONT00010	1
Set for presostat	PT	0100720005	1

## 1.8.TROUBLESHOOTING

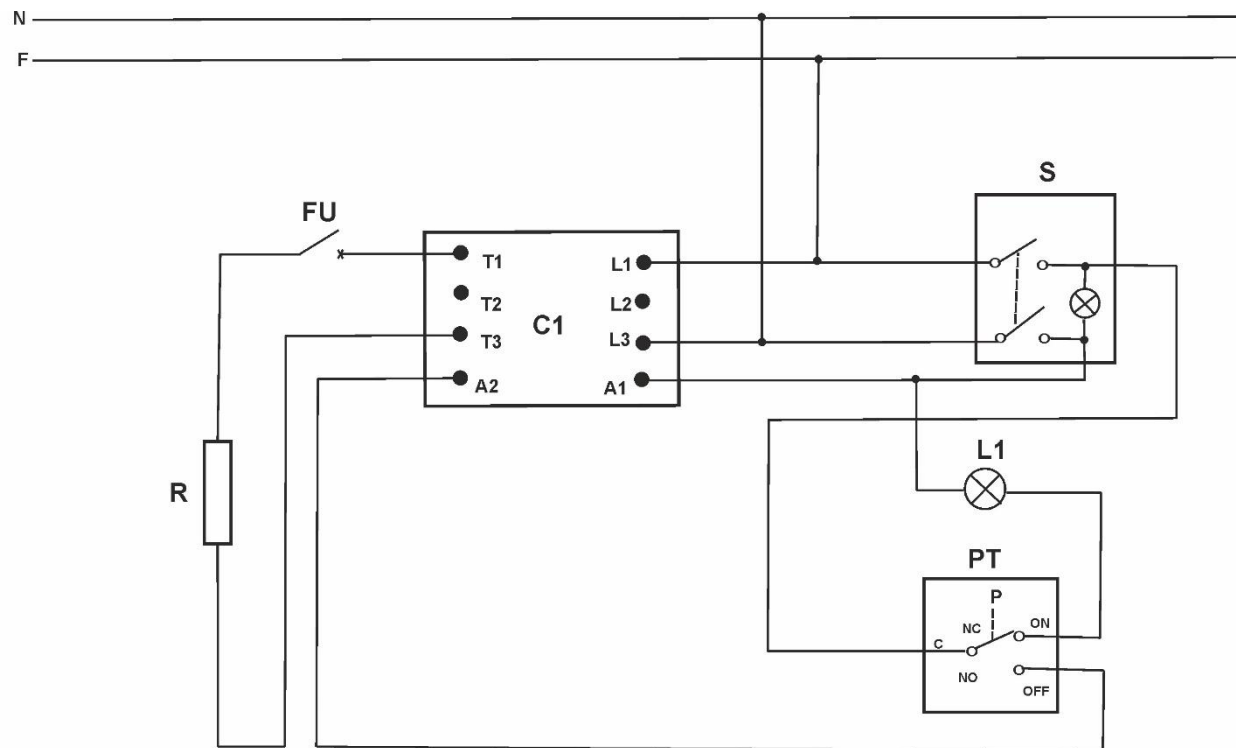PROBLEM	ACTION/S TO CARRY OUT
The green switch does not light when you switch on the water still	Check that the voltage of your electricity main is 230 V
The production of the water still is less than 4 L/H of distilled water	Check that the inlet water flow is 60 L/H Keep an appropriate level of water in the boiler Clean the heater
Low level of water in the boiler	Clean the lime quitting the obstruction in the hole between the level and boiler (please, see section 1.4, step 1)
The heater stops heating	Follow these steps by this order:  Check the state of the circuit breaker (FU).  The consumption of the heater is between 12.2 and 12.4 A  The yellow light must be off  The presostat must be commuting  The mini-contactor is working properly
The heater heats little	Clean the salt embedded on the heater
The heater is broken	Replace the heater

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## 1.10 WARRANTY

1. The product covered by this Certificate is protected for two (2) YEARS after the date of the purchase delivery note of the product.
2. During this period of time, in case of manufacturing fault of the product, the user has the right to its servicing, its replacement (except if one of this options was impossible or disproportionate), to have a reduction in the price and to the resolution of the contract (when the purchaser could not require the servicing or the replacement). The servicing will be free of charge for the user. All the servicing and replacements will be carried out in a reasonable period of time.
3. The breakdowns, flaws or caused damages in the product will be excluded as consequence of: abuses, improper use, exposure to adverse weather conditions, spilled batteries, blows, falls, bad or inadequate installation, power failure, waste by normal use, servicing carried out by different people indicated by us, as well as modification of this Certificate and / or another external causes to the product.
4. The glass parts of our distillers are excluded of this warranty.
5. This warranty will be automatically cancelled in case of wrong handling of the product or there is a breach of the manual or instructions regarding the use, assemble and operation.
6. The warranty does not affect to the rights of the customer according to the law.
7. In case the user needs technical help regarding the guaranteed product, the user will contact to the distributor.
8. With the exception of choice of law rules under Spanish private international law and the provisions of the UN Convention on Contracts for the International Sale of Goods (“UN-CISG”), the substantive law of Spain shall apply to all legal relationships and transactions established by this purchase agreement.
9. Please, should give copy of the delivery note or purchase invoice with the Warranty.